

PCR RIP B7

1- PCR mix

2,50 µl 10X buffer(Q BIOgene ; 1,5 mM MgCl₂)
 0,25 µl dNTP (20 mM total)
 0,15 µl primer B7a (30µM)
 0,15 µl primer B7b (30 µM)
 0,20 µl Taq polymerase (Q BIOgene ; 5U/µl)
 19,75 µl distilled water

2- PCR reaction :

23,0 µl mix
 2,0 µl DNA (200ng/µl)

3- PCR program :

5 min	94°C	
1 min	94°C	
1 min	60 °C	35 cycles
2 min	72°C	
5 min	72°C	

4- Primers:

- Primer B7a 5' GACACTGTTATACAGGGC 3'
 - Primer B7b 5' TGAAGCCATGGGCCACAC 3'

5- Amplification:

transgene : 850 bp